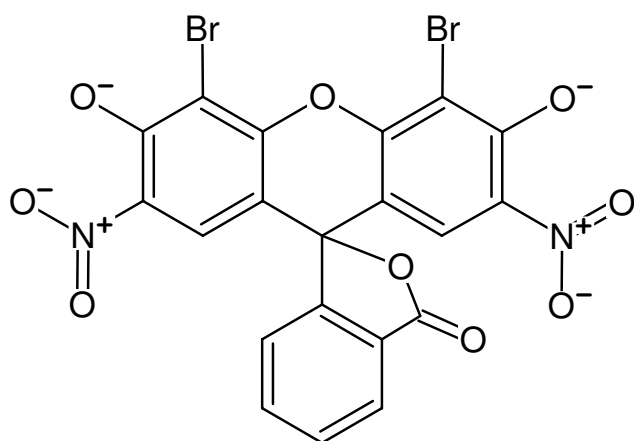



**JBS Bright Red**

Eosin Scarlet

Cat. No.	Amount
CO-304	300 µl

Na⁺Na⁺

Structural formula of JBS Bright Red

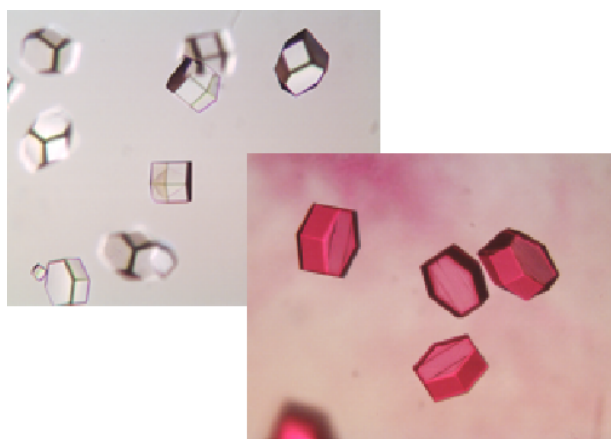

Unstained (A) and stained (B) protein crystals

For general laboratory use.**Shipping:** shipped at ambient temperature**Storage Conditions:** store at ambient temperature**Shelf Life:** 12 months**Molecular Formula:** C₂₀H₆Br₂N₂Na₂O₉**Molecular Weight:** 624.08 g/mol**CAS#:** 548-24-3**EC number:** 208-943-1**Concentration:** 1 mM**Applications:**

JBS Bright Red is a crystal dye used to stain macromolecular crystals, i.e. protein, peptide and nucleic acid crystals in order to differentiate them from small molecules and salt crystals.

Description:

Crystallization screening with high concentrations of precipitant and salt may lead to the formation of salt crystals. It is quite difficult to make a distinction between these false positives and true protein crystals.

Staining of crystals with appropriate dyes is a very straightforward method to differentiate between macromolecular crystals and salt crystals [1].

Protein and salt crystals differ substantially in their solvent content. Small crystal dyes, like JBS Bright Red, are able to permeate the solvent channels of a protein, thus coloring the protein red. In contrast, salt crystals are tightly packed and do not possess large solvent channels. They will therefore remain colourless.

Usage:

Simply add 0.5-1 µl of JBS Bright Red to the crystallization drop containing the crystals of interest.

Coloring Time:

JBS Bright Red colors protein crystals after a few minutes. Even if the color of the solution is only faintly red under the microscope, proteins will be stained within 5-15 min.

Very occasionally, it has been reported that protein crystals did not absorb crystal dyes [2].

Selected References:

- [1] Wilkosz *et al.* (1995) Preliminary characterization of EcoRI-DNA co-crystals: incomplete factorial design of oligonucleotide sequences. *Acta Cryst. D* **51**:938.
 [2] Eckert *et al.* (2003) Crystallization and preliminary X-ray analysis of



JBS Bright Red

Eosin Scarlet

Alicyclobacillus acidocaldarius endoglucanase CelA. *Acta Cryst. D* **59**:139.