

ReadiLink™ TCEP Removal Buffer

Ordering Information	Storage Conditions
Product Number: 5540 (1 mL, 1M)	-20°C

Standard Operating Protocol

1. Warm up ReadiLink™ TCEP removal buffer at room temperature before use.
2. Estimate the TCEP amount to be removed.
3. Add the ReadiLink™ TCEP removal buffer to a solution that contains residual TCEP at molar ratio of 100 : 1 TCEP Removal Buffer/TCEP (to be removed)
4. Incubate the solution (from Step 2) at room temperature from 30 to 60 minutes.
5. The solution is now ready to use for desired downstream assays.